


**Instruction:**

**Step 1.** Insert Adaptor 

**Step 2.** Connect HDMI cable from IMB08 to Monitor / TV, HDMI OUT 

**Step 3.** Connect HDMI cable from microscope to IMB08, HDMI IN 

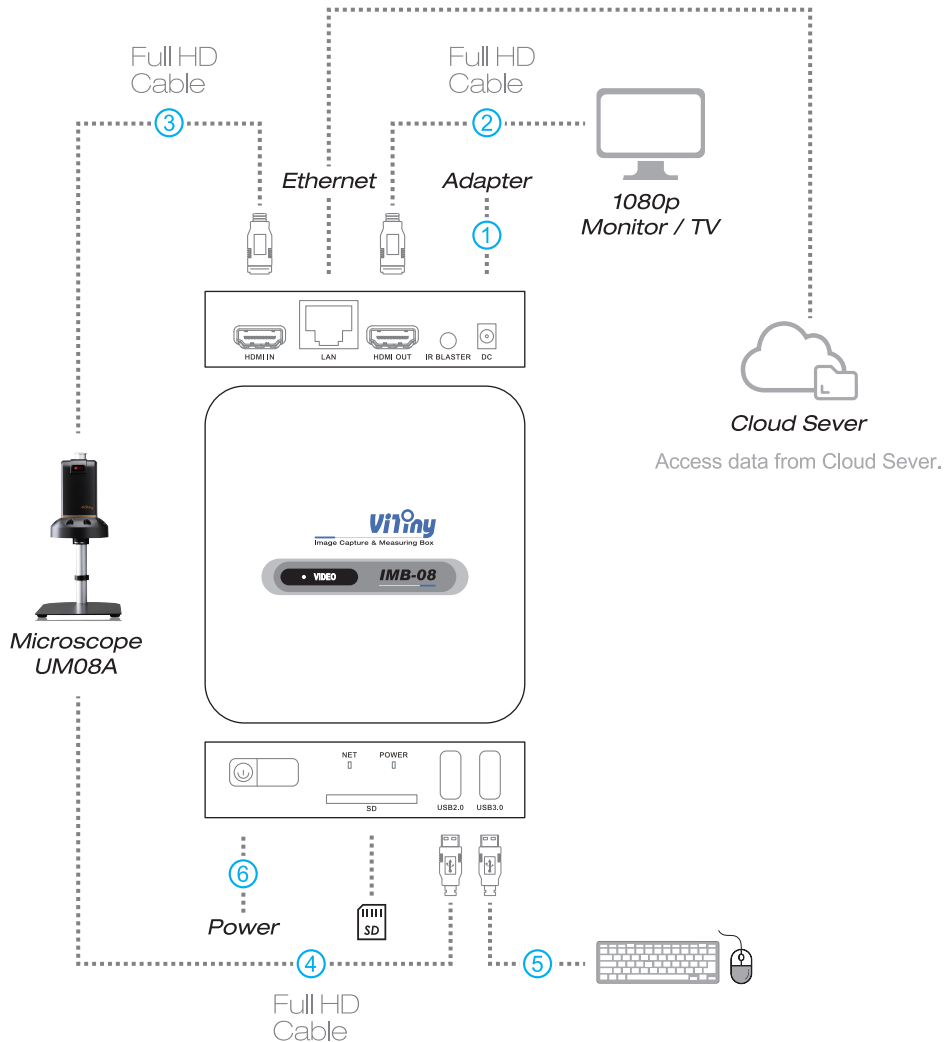

**Step 4.** Connect USB cable from microscope to IMB08 USB 2.0 

**Step 5.** USB 3.0 for external Mouse, Keyboard, USB Hub...etc. 

**Step 6.** Long press to power on. Short press for Restart and Shut Down mode. 


**Note :** SD Card slot : Store Image / Video.

# UmViewer App:

Long press to Power on




Click 「UmViewer」 App from Desktop

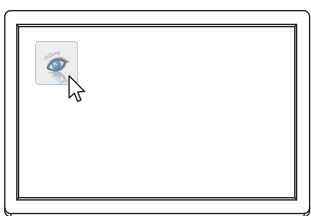

Function Tools

Preview Window

Picture / Video

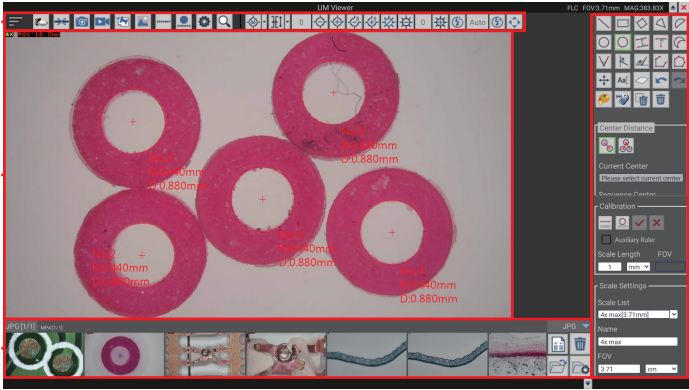

Measurement Tools

## Specification:

Internal Memory | 2GB DDR4

Internal Storage | 8GB

OS | Android 6.0

USB | USB2.0 x1  
| USB3.0 x1

HDMI | HDMI IN x1  
| HDMI OUT x1

DC Input | AC100 ~ 240V 50/60Hz

SD Slot | Max 32G

Video Output | HDMI

LAN | RJ45 10/100/1000M

Net Weight(g) | 212g

Size | 100x106x25(mm)

Power Adaptor | 5.0V-2A

## Package Contents:

| Item              | Q'ty |
|-------------------|------|
| Image Capture Box | 1    |
| HDMI Cable        | 1    |
| USB Cable         | 1    |
| Calibrator        | 1    |
| Adaptor           | 1    |